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User's Manual

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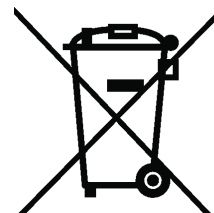
Contents

1	Introduction	5
1.1	Intended Use	5
1.2	Occupational Safety	5
2	Installing the Unit and Understanding its Functions	5
2.1	Installing and Checking out	5
2.2	Monitoring Unit Components	6
2.3	Understanding Operation	6
3	Operation	7
3.1	Normal Operation and Error Alert	7
3.2	Configuring the Settings	7
3.2.1	The Configuration Menu 1	8
3.2.2	Setting the Date and Time of Day	8
3.2.3	Selecting the Display Language	8
3.2.4	The Configuration Menu 2	8
3.2.5	Connecting a Cool Line Refrigerator (2.00.360C model version only)	8
3.2.6	Enabling an Audible Alarm in Case of Extraction Airflow Error	9
3.3	Potential-free ("Dry") Contacts	9
3.4	Setting the Pressure Difference	10
3.5	Operation on Battery, Switching on and off with an without Battery	11
3.6	Error and Menu Change Reporting	11
4	Technical Specifications	12
5	Troubleshooting	12
6	Contact Details	13

1 Introduction

Dear Customer:

Thank you very much for purchasing our product. We are confident that it will meet your expectations and perform satisfactorily in your daily routines.

The development, manufacture, and check of the ATEX-compliant exhaust airflow monitoring unit has been guided by our concern to offer a device characterized by superior operation safety and user-friendliness. The unit was manufactured and checked according to state-of-the-art technology and widely recognized safety regulations.

1.1 Intended Use

The exhaust air monitor has to be used in EX-free areas and is intended exclusively for the positive extraction of exhaust air from safety cabinets for storage installed indoors, zone 2 for gases belonging to temperature classes T1 - T4, gas type IIA or gas type IIB. The product is suitable only for fluid flows not transporting any particles or foreign bodies. The maximum temperature of extracted medium is 50°C (122°F).

Only such substances are acceptable for extraction that do not compromise the PP's connector sleeve. The relevant list of substances can be provided by the manufacturer in a digital format. Feel free to be in touch for any questions.

Caution

Any and all uses other than the above-described is prohibited and may cause hazardous situations, potentially danger of death, damage to the system and material property, and will cancel any and all warranty and liability for the product on the part of its manufacturer and distributor.

1.2 Occupational Safety

The installation and operation of the product must be performed in compliance with any and all nationally applicable rules and regulations for the storage of flammable liquids, the ISO 13857 standard, and any and all other applicable codes and standards. Modifications to the enclosure or the connector sleeve are prohibited.

Caution

The product must never be used if not totally installed, with the differential pressure meter properly fitted into the extraction system on site.

Electrical work on the product must only be performed when the exhaust air monitor was previously disconnected from power and/or locked against unwanted reapplication of power. This also holds for electrical work on the extraction system of which the monitor may be a part.

2 Installing the Unit and Understanding its Functions

2.1 Installing and Checking out

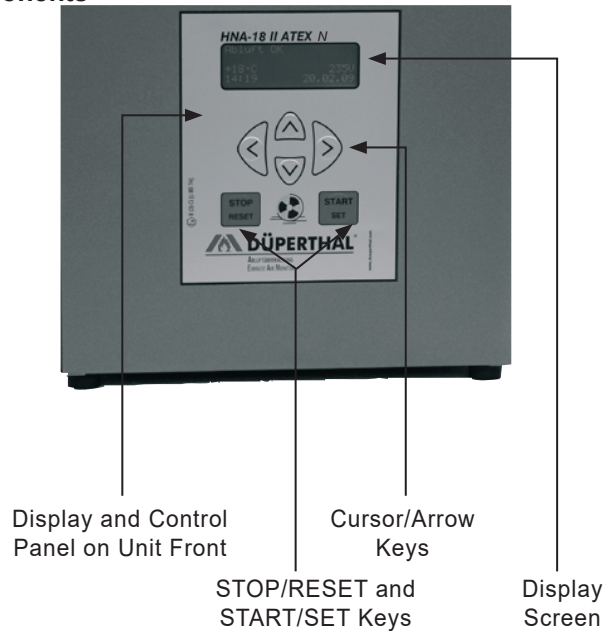
1. If needed or desired, connect the potential-free contacts (so-called "dry" contacts) to the terminal strip (see Sect. 3.3).
2. Connect the unit to the air extraction system on site:
Set the sensing sleeve with the differential pressure meter into the exhaust air line of the existing extractor and ensure that the con-

nection is tight. This is indispensable to ensure proper operation of the monitoring unit.

3. Connect the exhaust airflow monitor to power. Then wait a few moments for the monitor to initialize.
4. In some cases, it may be necessary to adapt the monitor to the existing extraction system. This involves mainly setting the differential pressure meter (see Sect. 3.4).

2.2 Monitoring Unit Components


Front View

i Where presses of the cursor keys are referenced below in this manual, they are symbolized by arrow characters:

- ← Key LEFT
- Key RIGHT
- ↑ Key UP
- ↓ Key DOWN

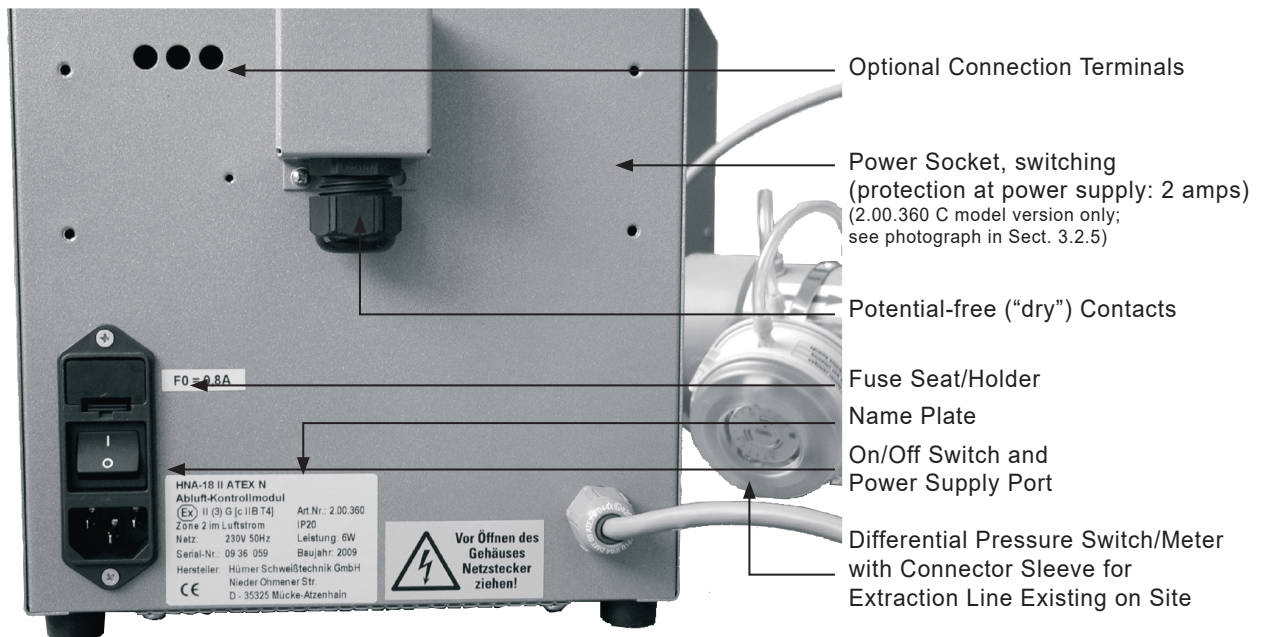Back View

2.3 Understanding Operation

After the On/Off switch was pressed to turn the module on, the airflow from the monitored system is constantly checked by the differential pressure meter inserted into the existing extraction line (see Figures in Sect. 2.2).

If an error condition develops for any reason, airflow through the line is restricted to a greater or lesser extent, and this causes the pressure difference measured in the suction hose to drop. If the pressure difference drops below the preset threshold, an audible alarm and a visual alarm are triggered.

If an error condition occurs, first the problem has to be cleared to re-enable proper airflow through the exhaust line. Then, as shown on the screen, the error and its correction have to be confirmed by pressing the START/SET key.

3 Operation

Caution

Be sure that all provisions of Sect. 1.1 and 1.2 above are strictly complied with for both the installation and the operation. Failure to do so may expose you to hazardous situations.

3.1 Normal Operation and Error Alert

When it works normally, the unit shows the so-called standard display (see Display 1) after it was switched on. The standard display shows the outside temperature, the detected power supply, and the date and time.

In case of an error condition, the reason for it is displayed. The example of Display 2 hints at a low voltage condition or power supply failure. An "Exhaust Air Error" may also be displayed. This error indicates another restriction, which is mechanical in most cases, of the flow of the fluid.

After an error was cleared, the unit indicates that the exhaust airflow has returned to normal. Still, the error and its successful correction have yet to be confirmed by pressing the START/SET key (see Display 3).

3.2 Configuring the Settings

Pressing the ⇐ and ⇒ keys simultaneously opens the configuration menus of the exhaust airflow monitor in which its parameters can be adjusted to your particular ways of operating the unit.

The 3 configuration menus are presented hierarchically and every level is accessible from the immediately lower level in the hierarchy. Whenever an operator wants to access one of the menu levels, either by using the above keypress combination or from the lower menu level, the airflow monitor requests entering an access code. The configuration options that are relevant for the operator are found on the two menu levels explained below (Configuration Menu 1 and Configuration Menu 2). Access to the hierarchically highest menu level, the so-called Supervisor Menu, is restricted to specifically trained technical staff.

When a menu is accessed, the access code applicable to the menu level in question is entered using the four arrow keys. After its input, the code has to be confirmed by pressing the START/SET key. The access code to access Configuration Menu 1 is **121**.

Use the ↑, ↓ arrow keys to select the desired configuration option. Use the ⇒ arrow key to toggle between "on" and "off" for that configuration option or to access a sub-menu for options where a symbolized arrow "→" is displayed (see Display 4).

Info

If a menu was accessed but no further action performed, the menu closes again automatically after approx. 1 minute, and no change is saved to memory.

If a menu was accessed and closed again, automatically or by the operator, it can be re-accessed without entering the access code again for 20 seconds after it was closed.

```
Exhaust Air OK
+23°C          231V
15:43          19.10.08
```

Display 1

```
Power Supply Error
+23°C          0V
15:59          19.10.08
```

Display 2

```
Exhaust Air OK
Confirm error!
+23°C          232V
16:06          19.10.08
```

Display 3

```
* ENTER MENU CODE *
***
```

Display 4a

```
** Config Menu 1 **
Config Menu 2    ->
>Date/Time       ->
Language         ->
```

Display 4b

3.2.1 The Configuration Menu 1

At the level of Configuration Menu 1 (see Display 4b), the following settings can be configured after the access code **121** was entered.

“Configuration Menu 2 →” means that by pressing the ⇨ key, the user can access the level of Configuration Menu 2, provided the access code corresponding to that menu level is entered.

“Date/Time →” means that by pressing the ⇨ key, the user can access a sub-menu for setting the date and the time of day.

“Language →” means that by pressing the ⇨ key, the user can access a sub-menu for selecting the display language.

When the Configuration Menu 1 was accessed, the currently displayed settings are saved to memory by pressing the START/SET key.

3.2.2 Setting the Date and Time of Day

When the “Date/Time” sub-menu was selected, the screen changes to the display reproduced in Display 5.

The date and the time of day can be set using the cursor keys. Move the cursor (underline symbol) to the digit you want to change by pressing the ⇐ and ⇒ keys, change it by pressing the ↑ and ↓ keys. The set date and the set time must not be in the past.

Press the START/SET key to confirm and save your settings.

```
*** Date/Time ***
19.10.08      17:01
```

Display 5

3.2.3 Selecting the Display Language

When the “Language” sub-menu was selected, the screen changes to the display reproduced in Display 6, where the arrow symbol indicates the currently selected language. Use the arrow keys ↑ and ↓ to select the desired language for the display screen, and confirm your selection by pressing the START/SET key.

```
Deutsch
>English
Français
```

Display 6

3.2.4 The Configuration Menu 2

At the level of Configuration Menu 2 (see Display 7b), which is accessible from Configuration Menu 1, the following settings can be configured. The access code required is **41531**. It has to be entered when access to this menu is requested.

“Supervisor Menu →” means that by pressing the ⇨ key, the user can access the level of Supervisor Menu, which contains system parameters that are unimportant for standard operation. Please contact the manufacturer for further information on this matter.

“Exhaust Monitor on” means that both power supply and airflow are monitored, “off” that only the power supply is monitored, no monitoring of extraction flow takes place. Accordingly, when exhaust airflow monitoring is off, only power supply errors will be detected.

“Cool Line on” means that the power socket at the back of the enclosure, which can power a Cool Line refrigerator cabinet, is switched on, “off” that the socket is dead and, hence, power to the refrigerator is switched off. (The power socket and this menu option are available only on units of the 2.00.360C model version; see Sect. 3.2.5.)

“Airflow Signal →” means that pressing the ⇨ key opens a sub-menu for enabling and setting an audible alarm to sound with the error message in the event of an exhaust airflow error (see Sect. 3.2.6).

When the Configuration Menu 2 was accessed, the currently displayed settings are saved to memory by pressing the START/SET key.

```
* ENTER MENU CODE *
*****
```

Display 7a

```
** Config Menu 2 **
Supervisor Menu ->
Exhaust Monitor on
>Cool Line      on
```

Display 7b

3.2.5 Connecting a Cool Line Refrigerator (2.00.360C model version only)

The custom-designed power socket at the back of the enclosure (see the

photograph on the next page), which is designed for connecting a Cool Line series refrigerator to the unit, can be enabled or disabled with the “Cool Line” option of Configuration Menu 2. When it is disabled, the refrigerator is switched off and there is a “Cool Line disabled” message on the standard display, alternating with the status message “Exhaust Air OK” on the first line. Acceptable specifications of a refrigerator are 230 V AC, 50 Hz, max. 5 A.

Important

The 2.00.360 C model version of the exhaust air control unit ships with a dedicated adapter cable matched to the custom-designed power socket (see the photographs to the right). Power may be supplied to the cabinet **exclusively** with the factory-installed socket and shipped, dedicated cable.

Important

As soon as an error condition is detected, power to the power socket is switched off, which is equivalent to turning the refrigerator off. The same holds when the exhaust air monitoring is disabled in the Configuration Menu 2.

Info

When enabled, the special power socket at the back of the exhaust air control unit checks whether it receives a loop-back signal from the connected storage cabinet. If it is enabled, but does not detect a loop-back signal, a message “Error Cool Line” is displayed at the seconde line of the screen.

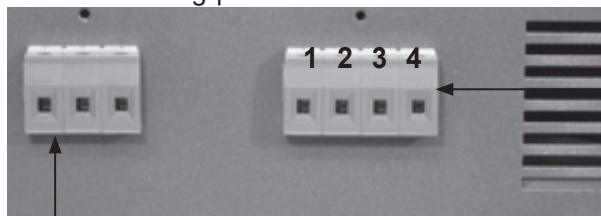
3.2.6 Enabling an Audible Alarm in Case of Extraction Airflow Error

When the “Airflow Signal” submenu was selected, the ↑, ↓ cursor keys are used to select the desired behavior of the audible alarm signal that sounds with an extraction airflow error indicated on the screen. The selected item is saved and applied when you press START/SET.



The “No Delay” item means that the alarm will sound immediately when an error message appears on the screen. By selecting “Off,” you disable the alarm sound altogether. The time items (30 sec, 1 min, 1.5 min, 2 min, 3 min), when selected, will trigger the sound with a delay of 30 sec or 1 min etc. from the moment the error message is displayed.

3.3 Potential-free (“Dry”) Contacts

An error condition causes two normally open, potential-free contacts to be closed; for individually locating errors, their closure can be recorded at the switching panel from the backside of the enclosure (see above).

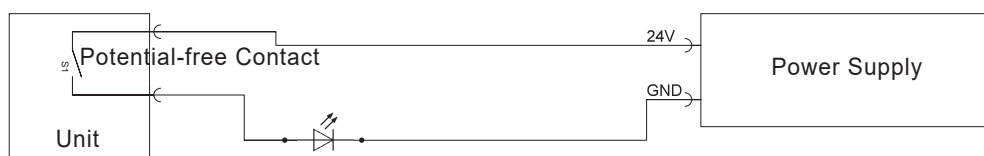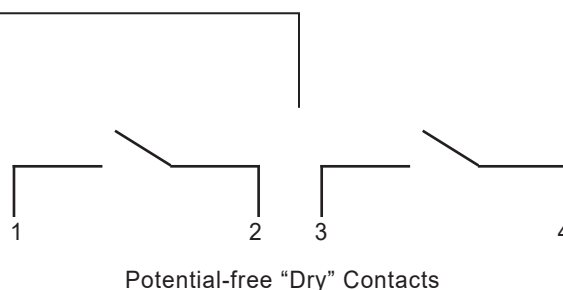Prepared for the Installation of a Mechanical or Inductive Position Switch

Confirming the error and its correction by pressing the START/SET key (see Sect. 3.1) causes the two contacts to be reset (i.e., opened).

Important

The potential-free circuit (only one contact reproduced in the diagram) is designed for **max. 24 V, 2 A**. If these values were exceeded, the unit may be irrevocably destroyed. Additionally, any such damage would not be covered by the product warranty.

3.4 Setting the Pressure Difference

At a metering orifice, the suction line from the storage cabinet is equipped with a pitot probe. The pressure difference it senses is measured by a differential pressure meter via the two transparent tubes routed to it. It rises as the airflow through the suction-side extraction line on site increases. With restricted or blocked airflow, the pressure difference sensed by the probe drops. If it drops below the threshold set at the differential pressure meter, a signal that an error condition has developed is sent into the enclosure for further processing via the cable in the blue protective conduit. For technical reasons (hysteresis), the threshold from which the monitor cancels the error condition when the pressure difference rises again, is a little higher than the threshold that triggers an error. Therefore, make sure at all times that the settings for the pressure difference and the fan speed of the air extraction system on site are matched to each other in such a way that both an error is indicated when a restriction is detected, and the error is canceled by the unit when the restriction is cleared. There is no upper threshold for a maximum pressure difference.

The value of the pressure difference threshold can be set at the meter from 20 Pa to 300 Pa. This is done by simply removing the transparent cover and using a screwdriver in the provided slot to turn the selector. The factory-set default threshold of the pressure difference is 70 Pa.

Setting the pressure difference threshold for monitoring the airflow requires the following steps:

1. Connect the exhaust air monitoring unit properly to the extraction system (outlet nipple) on site. Pay attention to the installation of the sleeve of the differential pressure meter; it has to tee into the line so the airflow passes through it in the direction of the arrow.
2. Set the fan speed to the desired extraction airflow (see the operator instructions of the fan system on site).

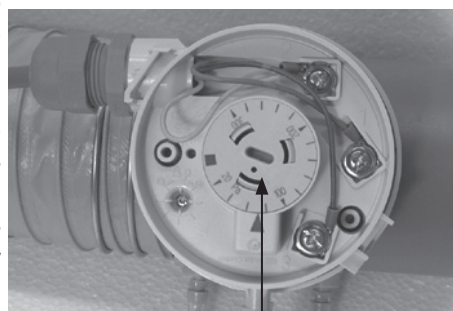
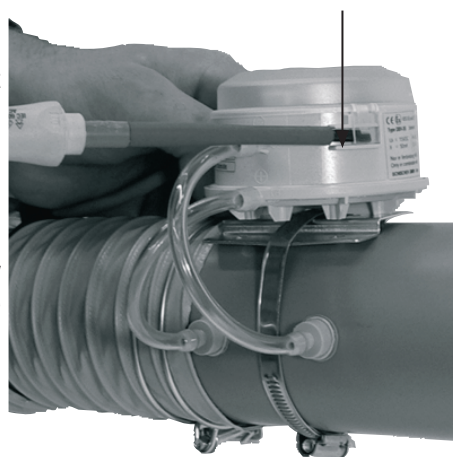
Caution

In a cabinet for flammable liquids, the extraction airflow has to be sufficient to replace the interior volume of the cabinet 10 times every hour.

3. Loosen the plastic cover of the differential pressure meter on the suction hose (refer to the photograph to the right).
4. Using an appropriate screwdriver in the turning selector at the center of the differential pressure meter, set the threshold to make "Exhaust OK" appear on the screen (reduce pressure difference). If no error is displayed on the screen at the start, the pressure difference setting may first be increased until an error is displayed, then slowly reduced again. The higher the pressure difference threshold is set, the quicker / more sensitive the meter will respond by throwing an error message.
5. Using appropriate means on site (e.g. by covering the exhaust outlet with your hand), interrupt the extraction airflow temporarily.
6. If set correctly, the unit will trigger the audible alarm and display the error message on the screen. After the airflow interruption

Latch clips to hold/remove the plastic cover

Selector for setting the pressure difference threshold

is cleared, the audible alarm should stop and the "Confirm error" message should appear on the screen.

7. Reinstall the plastic cover on the differential pressure meter.
8. Confirm the error message on the screen by pressing the START/SET key to re-open the dry contacts.

3.5 Operation on Battery, Switching on and off with an without Battery

In the event of power supply failure, the control unit can be operated on the internal battery (accumulator) for approx. 3 hours.

This is why, provided the battery is sufficiently charged, pressing the On/Off switch will not turn the unit off. When the switch was pressed, the battery will continue to supply power to it. To turn the monitoring unit off definitely, i.e. to disconnect power to it from both the mains network and the battery, hold down the STOP/RESET key for approx. 5 seconds when the standard display shows on the screen. This causes the message "Please disconnect power to unit!" to appear. Do as you are asked to: switch the unit off using the On/Off switch and unplug the mains plug. Wait a short time for the unit to switch itself off.

If the unit is switched off without holding the STOP/RESET key down first, this counts as a power supply failure. It will trigger an error message and close the dry contacts (see Sect. 3.1 and 3.3). In case the On/Off switch was pressed first to switch the unit off, the unit will indeed switch itself off when the STOP/RESET key is then held down. However, the error message that the switch caused has to be confirmed when it is switched on again.

If the battery is charged to a sufficient level, it is possible to switch the unit on while it is not connected to mains power, by pressing the START/SET key. If the battery voltage drops below a factory-set threshold, the monitoring unit switches itself off automatically. Turning on and running the unit without mains power will trigger a power supply error message as well, which will have to be confirmed.

3.6 Error and Menu Change Reporting

Error conditions and changes applied to the configuration menus are recorded and saved to memory. They can be viewed after pressing the ⇐ arrow key in the standard display (see Display 8). You can scroll through the displayed listing by pressing the ↑ and ↓ arrow keys. The maximum number of entries saved to memory and displayed in this listing in chronological order is 50. When the maximum number is reached, any further report will overwrite the oldest one in memory. Deleting the entries in memory is possible from the Supervisor Menu.

A report entry contains information on the type of entry, i.e. error or menu change, the date, and the time of day. The title line of the listing shows the number of the report currently looked at (before the slash) and the total number of entries in memory (after the slash). If the entry is an error report, the error code (E1 or E2) is shown. These error codes have the following meaning:

- E1 Power Supply Error
- E2 Exhaust Air Error

```
*   Report 018/42   *
E1 22.06.08 13:32
>M 22.06.08 13:41->
E2 24.06.08 08:09
```

Display 8

The letter "M" stands for the report of a modification of the configuration menus. As is indicated by the symbolized arrow to the right, pressing the ⇨ arrow key, when this entry is selected, will show additional information. This information is a list of all settings of the configuration menus. All the options are given with the indication of their status, changed or

unchanged at the moment the modification record was generated. The record for a modification is generated only when the operator quits the configuration menus, and it includes all settings, whether changed at that point in time or remained as is.

The following overview is meant to assist you in interpreting the change reports.

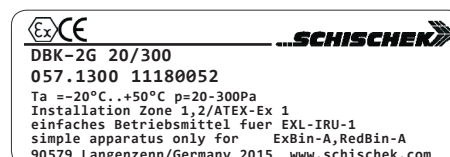
Date/Time <> Date/time before do not equal after change
 Language = Set language before equals after change
 Exh Monitor 1->0 The exhaust monitoring setting was set from "on" (1) to "off" (2)
 Cool Line 1->1 The setting for the refrigerator socket was left at "on" (with 2.00.360C models only)

The report screen is quit by pressing the STOP/RESET key, or it returns automatically to the standard display after a short time of inactivity.

4 Technical Specifications

Power Supply Voltage 230 V
 Frequency 50 Hz
 Current Rating 0.03 A
 Weight approx. 3.5 kg
 Dimensions (enclosure edges) 215 x 175 x 220 mm
 Ingress Protection IP20
 ATEX Rating EX II (3) G [c II B T4]

see also the component name plate, as to the right

5 Troubleshooting

In the course of using the appliance, malfunctions may occur. The table below is meant to help troubleshoot and correct malfunctions. If a malfunction cannot be traced to one of the possible causes or the correction given does not solve the issue, then use of the appliance has to be discontinued and it has to be returned for inspection and repair.

Malfunction	Possible Cause	Verification of Cause	Correction of Malfunction
Appliance does not run	On/Off Switch Power Supply Cord Overvoltage Switch Power Socket at the Back (Cool-Line cabinet)	- Is the appliance switched on at the On/Off switch? - Is the power supply cord damaged? - Is the power supply cord properly connected? - Are the fuses inside the appliance o.k.? - Is a device connected that may consume current at more than 5 A?	- Switch the appliance on at the On/Off switch. - Replace the supply cord with a new one. - Connect the supply cord to the intended socket. - Replace the fuses (2 x 0.8 A slow blow). - Disconnect this device from the socket and check the fuses and replace them as needed (2 x 5 A slow blow).
Display screen does not work	Display Electronics/ Electronics	Malfunction cannot be corrected by user.	Return the appliance to the manufacturer/ seller for inspection/repair.
Power Supply Error (E1) shows on screen	Possible Cause, Verification, and Correction as under "Appliance does not run" above.		

Malfunction	Possible Cause	Verification of Cause	Correction of Malfunction
Exhaust Airflow Error (E2) shows on screen	Differential Pressure Meter/Switch	- Are the measuring tubes connecting the sleeve to the meter/switch properly connected? - Are the measuring tubes connecting the sleeve to the meter/switch damaged? - Are the measuring tubes connecting the sleeve to the meter/switch connected to the correct nipples? - Is the differential pressure meter/switch properly set? - Is the extraction line on site connected tightly to the sleeve? - Is the differential pressure meter/switch wired properly?	- Connect the measuring and connecting tubes to both the meter/switch and the sleeve in the extraction line. - Replace the measuring and connecting tubes with new ones. - Connect the tubes as intended: nipple closer to the cabinet of the sleeve to P1 (+) of the meter/switch, nipple closer to the appliance of the sleeve to P2 (–) of the meter/switch. - Open the differential pressure meter/switch and set it properly (see Sect. 3.4). - Tightly set the sleeve into the extraction line on site, for instance using the clamps at both sleeve ends. - Open the differential pressure meter/switch and wire it correctly: blue wire to 1, yellow to 2, red to 3. - Return the appliance to the manufacturer/seller for inspection/repair.
Despite irregular extraction airflow no airflow error (E2) on screen	Switching Amplifier	Is a clapping noise heard when the differential pressure switch cuts in or out?	
	Configuration Menu 2 Differ. Pressure Meter/Switch	- Is the feature for monitoring the extraction airflow turned off? - Is the differential pressure meter/switch properly set?	- Set the menu option for the exhaust airflow monitor to “on”. - Open the differential pressure meter/switch and set it properly (see Sect. 3.4).

6 Contact Details

Düperthal Sicherheitstechnik GmbH & Co. KG
 Frankenstr. 3 Tel.: +49 6188 9139 0
 63791 Karlstein, Germany Fax: +49 6188 9139 121
 Web: www.dueperthal.com E-mail: info@dueperthal.com

We reserve the right to change technical specs of the unit without prior notice.

Pursuant to the directive 2012/19/EU on Waste Electrical and Electronic Equipment (so-called WEEE Directive), equipment that was manufactured or distributed by ourselves may be returned to us. To discuss the exact procedure we will follow, please contact us with the details above.

We also declare that the equipment manufacture conforms to the directive 2011/65/EU on the restriction of the use of certain hazardous substances in electrical and electronic equipment (so-called RoHS Directive).

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