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## User's Manual

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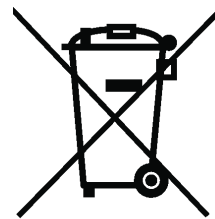
# Contents

1	Introduction .....	5
1.1	Explaining Icons Used.....	5
1.2	Intended Use.....	5
1.3	Occupational Safety .....	5
2	Installing the Unit and Understanding its Functions .....	6
2.1	Installing and Checking out .....	6
2.2	Monitoring Unit Components.....	6
2.3	Understanding Operation .....	7
3	Operation .....	7
3.1	Normal Operation and Error Alert.....	7
3.2	Configuring the Settings.....	8
3.2.1	The Configuration Menu 1.....	8
3.2.2	Setting the Date and Time of Day .....	8
3.2.3	Selecting the Display Language.....	9
3.2.4	The Configuration Menu 2.....	9
3.2.5	Setting the Speed of the Fan .....	9
3.2.6	Connecting a Cool Line Refrigerator (2.00.321 C model version only) .....	10
3.2.7	Enabling an Audible Alarm in Case of Extraction Airflow Error .....	10
3.3	Potential-free (“Dry”) Contacts .....	11
3.4	Setting the Pressure Difference .....	11
3.5	Operation on Battery, Switching on and off with and without Battery.....	12
3.6	Online Monitoring over the Internet with DÜPERTHAL connect .....	13
3.7	Error and Menu Change Reporting .....	13
4	Technical Specifications .....	14
5	Troubleshooting.....	14
6	Contact Details.....	15

## 1 Introduction

Dear Customer:

Thank you very much for purchasing our product. We are confident that it will meet your expectations and perform satisfactorily in your daily routines.

The development, manufacture, and check of the ATEX-compliant exhaust airflow monitoring unit with fan has been guided by our concern to offer a device characterized by superior operation safety and user-friendliness. The unit was manufactured and checked according to state-of-the-art technology and widely recognized safety regulations.

### 1.1 Explaining Icons Used

The following expressions and icons are used in this User's Manual to refer to safety-related issues:

Caution

This icon indicates that non-compliance may result in a hazardous situation that possibly causes bodily injury or material damage.

Important

This icon indicates important messages related to the correct use of the product. Non-compliance may cause problems of operation and damage to the product.

Info

This icon indicates tips and useful information for using the product more efficiently and more economically.

### 1.2 Intended Use

The exhaust air monitor with fan has to be used in EX-free areas and is intended exclusively for the positive extraction of exhaust air from safety cabinets for storage installed indoors, zone 2 for gases belonging to temperature classes T1 - T4, gas type IIA or gas type IIB. The product is suitable only for fluid flows not transporting any particles or foreign bodies. The maximum temperature of extracted medium is 50°C (122°F).

Only such substances are acceptable for extraction that do not compromise the PPS connector sleeve, the PVC suction hose, and the fan. The relevant list of substances can be provided by the manufacturer in a digital format. Feel free to be in touch for any questions.

Caution

Any and all uses other than the above-described is prohibited and may cause hazardous situations, potentially danger of death, damage to the system and material property, and will cancel any and all warranty and liability for the product on the part of its manufacturer and distributor.

From the model version capable of radio communication (P/N. 2.00.322), a wireless connection can be established to DÜPERTHAL connect. This requires a Bridge to be present as a gateway in the radio-covered area.

### 1.3 Occupational Safety

The installation and operation of the product must be performed in compliance with any and all nationally applicable rules and regulations for the storage of flammable liquids, the ISO 13857 standard, and any and all other applicable codes and standards. The design of the product is such that the fan blades are off-touch. Modifications to the enclosure or the suction and exhaust hoses that may make the fan blades accessible to the touch when the unit is in normal use are prohibited.

Caution

The product must never be used if not totally installed, with properly set suction and exhaust hoses. Electrical work on the product must only be performed when the exhaust air monitor was previously disconnected from power and/or locked against unwanted reapplication of power. This also holds for electrical work on the extraction system of which the monitor may be a part.

## 2 Installing the Unit and Understanding its Functions

### 2.1 Installing and Checking out

1. If needed or desired, connect the potential-free contacts (so-called “dry” contacts) to the terminal strip (see Sect. 3.3).
2. Connect the unit to the air extraction system on site:  
 Connect the existing ventilation system to the unit, e.g., with a flexible sleeve fitting, making sure no strain is imposed on the connection, and make a tight connection using an appropriate hose clamp. This is indispensable to ensure proper operation of the extraction fan. As a general rule, be sure that the system is set up with air lines that are as short as possible.

Caution

In no circumstances, it is acceptable to transfer the air extracted by the blower fan back into the work room.

3. Connect the exhaust airflow monitor to power. Then wait a few moments for the monitor to initialize.
4. In some cases, it may be necessary to adapt the monitor to the existing extraction system. This involves mainly setting the differential pressure meter and/or the speed of the fan (see Sect. 3.4).

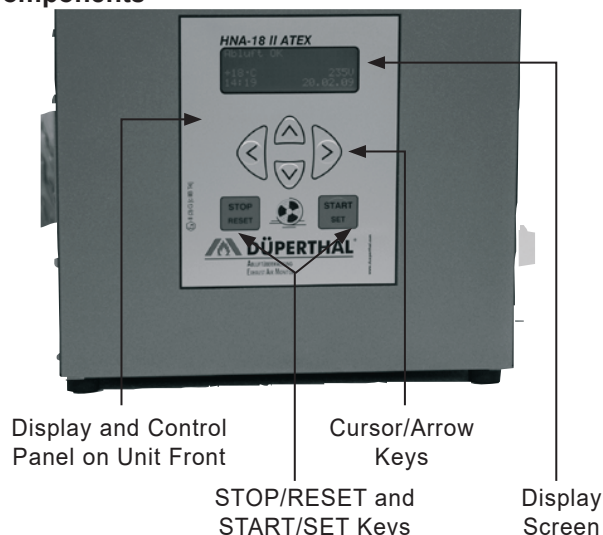
Info

Be aware that the actually existing pressure difference is directly dependent upon the speed with which the fan turns. Therefore, if you are authorized to change the fan speed, bear in mind that the fan speed and the pressure difference threshold have to be matched to each other. If this were not the case, there would be a risk of getting false alarms.

5. Register, if you so desire, the unit with the DÜPERTHAL connect system for monitoring the operation online over the internet (possible only with the model version capable of radio communication, P/N. 2.00.322; see Sect. 3.6).

### 2.2 Monitoring Unit Components


Front View

Where presses of the cursor keys are referenced below in this manual, they are symbolized by arrow characters:

- ← Key LEFT
- Key RIGHT
- ↑ Key UP
- ↓ Key DOWN

## Back View
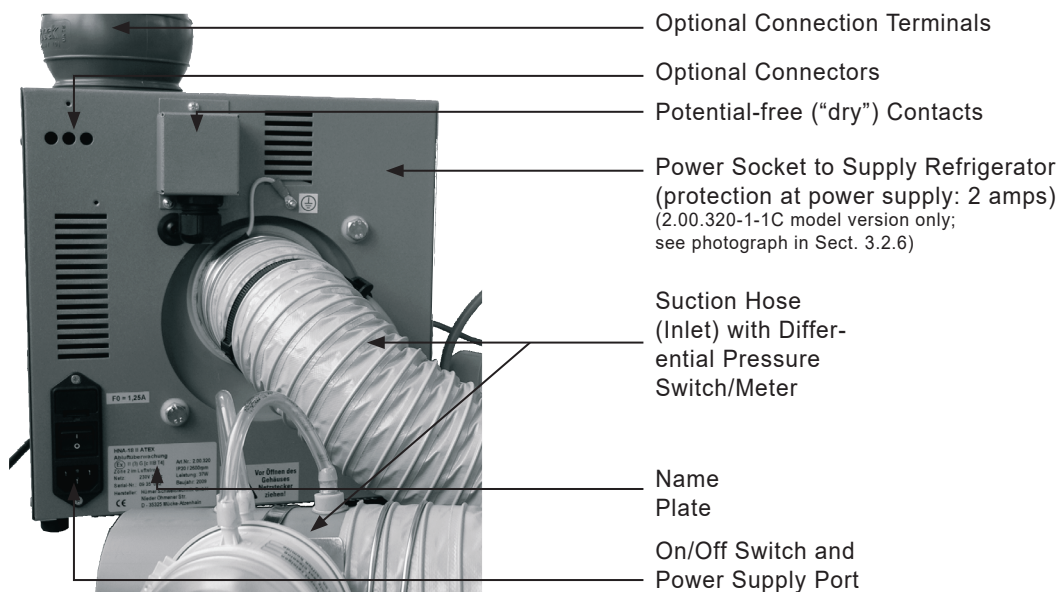
## 2.3 Understanding Operation

After the On/Off switch was pressed to turn the unit on, the fan located inside the enclosure of the air flow monitor draws air from the monitored system through the suction hose connected to the inlet nipple and extracts the air from the system through the blower nipple on the enclosure (see Figures in Sect. 2.2).

If an error condition develops, airflow through the line is restricted to a greater or lesser extent, and this causes the pressure difference measured in the suction hose to drop. If the pressure difference drops below the set threshold, an audible alarm and a visual alarm are triggered.

If an error condition occurs, first the problem has to be cleared to re-enable proper airflow through the exhaust line. Then, as shown on the screen, the error and its correction have to be confirmed by pressing the START/SET key.

## 3 Operation

Caution

Be sure that all provisions of Sect. 1.2 and 1.3 above are strictly complied with for both the installation and the operation. Failure to do so may expose you to hazardous situations.

### 3.1 Normal Operation and Error Alert

When it works normally, the unit shows the so-called standard display (see Display 1) after it was switched on. The standard display shows the outside temperature, the detected power supply, and the date and time.

```
Exhaust Air OK
+23°C                231V
15:43                19.10.08
```

Display 1

In case of an error condition, the reason for it is displayed. The example of Display 2 hints at a low voltage condition or power supply failure. An "Exhaust Air Error" may also be displayed. This error indicates another restriction, which is mechanical in most cases, of the flow of the fluid.

```
Power Supply Error
+23°C                0V
15:59                19.10.08
```

Display 2

After an error was cleared, the unit indicates that the exhaust airflow has returned to normal. Still, the error and its successful correction have yet to be confirmed by pressing the START/SET key (see Display 3).

```
Exhaust Air OK
Confirm error!
+23°C                232V
16:06                19.10.08
```

Display 3

Important

The fan has to be checked at regular intervals. If it is dirty, damaged or shows signs of corrosion, it has to be cleaned or

the defects have to be corrected, if needed by returning the blower box for repair to the seller or manufacturer.

### 3.2 Configuring the Settings

Pressing the ⇐ and ⇒ keys simultaneously opens the configuration menus of the exhaust airflow monitor in which its parameters can be adjusted to your particular ways of operating the unit.

The 3 configuration menus are presented hierarchically and every level is accessible from the immediately lower level in the hierarchy. Whenever an operator wants to access one of the menu levels, either by using the above keypress combination or from the lower menu level, the airflow monitor requests entering an access code. The configuration options that are relevant for the operator are found on the two menu levels explained below (Configuration Menu 1 and Configuration Menu 2). Access to the hierarchically highest menu level, the so-called Supervisor Menu, is restricted to specifically trained technical staff.

When a menu is accessed, the access code applicable to the menu level in question is entered using the four arrow keys. After its input, the code has to be confirmed by pressing the START/SET key. The access code to access Configuration Menu 1 is **121**.

Use the ↑, ↓ arrow keys to select the desired configuration option. Use the ⇒ arrow key to toggle between “on” and “off” for that configuration option or to access a sub-menu for options where a symbolized arrow “->” is displayed (see Display 4).

If a menu was accessed but no further action performed, the menu closes again automatically after approx. 1 minute, and no change is saved to memory.

If a menu was accessed and closed again, automatically or by the operator, it can be re-accessed without entering the access code again for 20 seconds after it was closed.

#### 3.2.1 The Configuration Menu 1

At the level of Configuration Menu 1 (see Displays 4, 5), the following settings can be configured after the access code **121** was entered.

“Configuration Menu 2 -> ” means that by pressing the ⇒ key, the user can access the level of Configuration Menu 2, provided the access code corresponding to that menu level is entered.

“Date/Time -> ” means that by pressing the ⇒ key, the user can access a sub-menu for setting the date and the time of day.

“Language -> ” means that by pressing the ⇒ key, the user can access a sub-menu for selecting the display language.

When the Configuration Menu 1 was accessed, the currently displayed settings are saved to memory by pressing the START/SET key.

#### 3.2.2 Setting the Date and Time of Day


When the “Date/Time” sub-menu was selected, the screen changes to the display reproduced in Display 6.

The date and the time of day can be set using the cursor keys. Move the cursor (underline symbol) to the digit you want to change by pressing the ⇐ and ⇒ keys, change it by pressing the ↑ and ↓ keys. The set date and the set time must not be in the past.


Press the START/SET key to confirm and save your settings.

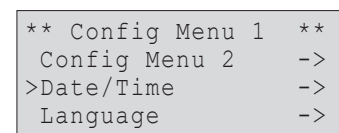Display 4


Display 5

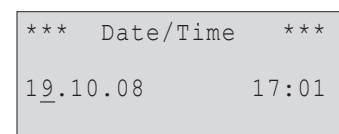Display 6

### 3.2.3 Selecting the Display Language

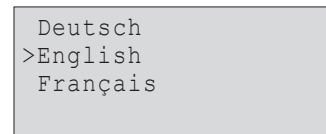When the “Language” sub-menu was selected, the screen changes to the display reproduced in Display 7, where the arrow symbol indicates the currently selected language. Use the arrow keys  $\uparrow$  and  $\downarrow$  to select the desired language for the display screen, and confirm your selection by pressing the START/SET key.

Display 7

### 3.2.4 The Configuration Menu 2

At the level of Configuration Menu 2 (see Displays 8, 9), which is accessible from Configuration Menu 1, the following settings can be configured. The access code required is **41531**. It has to be entered when access to this menu is requested.

“Supervisor Menu  $\rightarrow$ ” means that by pressing the  $\Rightarrow$  key, the user can access the level of Supervisor Menu, which contains system parameters that are unimportant for standard operation. Please contact the manufacturer for further information on this matter.

“Exhaust Monitor on” means that both power supply and airflow are monitored, “off” that only the power supply is monitored, no monitoring of extraction flow takes place. Accordingly, when exhaust airflow monitoring is off, only power supply errors will be detected.


“Fan Speed  $\rightarrow$ ” means that by pressing the  $\Rightarrow$  key, the user can access a sub-menu for setting the speed at which the fan should turn.

“Cool Line on” means that the power socket at the back of the enclosure, which can power a Cool Line refrigerator cabinet, is switched on, “off” that the socket is dead and, hence, power to the refrigerator is switched off. (The power socket and this menu option are available only on units of the 2.00.321 C model version; see Sect. 3.2.6.)

“Airflow Signal  $\rightarrow$ ” means that pressing the  $\Rightarrow$  key opens a sub-menu for enabling and setting an audible alarm to sound with the error message in the event of an exhaust airflow error (see Sect. 3.2.7).

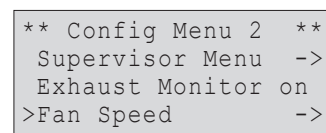Display 8


Display 9

When the Configuration Menu 2 was accessed, the currently displayed settings are saved to memory by pressing the START/SET key.

### 3.2.5 Setting the Speed of the Fan

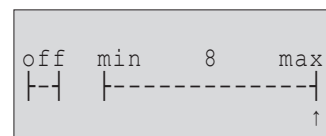When the “Fan Speed” sub-menu was selected, the screen changes to the display reproduced in Display 10. The currently set speed of the fan is indicated by the cursor arrow sign. When it is shipped ex works, the unit is set to maximum fan speed.

Press the  $\Leftarrow$  or  $\Rightarrow$  key to move the cursor to the left or the right, thereby reducing or increasing the fan speed. To turn the fan off, move the cursor once more to the left, pressing the  $\Leftarrow$  key, when it is on the minimum position. To set the new speed, press the START/SET key.

Display 10

Important

For increased safety, the STOP/RESET key is disabled while a modification of the fan speed setting is in progress, in order to avoid any inadvertent change. You will have to positively confirm the currently displayed value by pressing START/SET, to be able to quit this input screen.


Info

The information on fluid velocity and volume flow rate in the table below relate to the following setup:

- Inside diameter of measuring tube 70.5 mm
- Actual measured length approx. 3 m

Level	0	1	2	3	4	5	6	7	8	9	10
Fluid Velocity, m/s	0	1.42	1.65	1.83	2.06	2.32	2.50	2.74	2.98	3.15	3.41
Volum. Flow Rate, m³/h	0	20.0	23.2	25.7	29.0	32.6	35.0	38.5	41.9	44.3	47.9

Level	11	12	13	14	15
Fluid Velocity, m/s	3.48	3.85	4.07	4.34	4.51
Volum. Flow Rate, m³/h	48.9	54.0	57.2	60.9	63.5

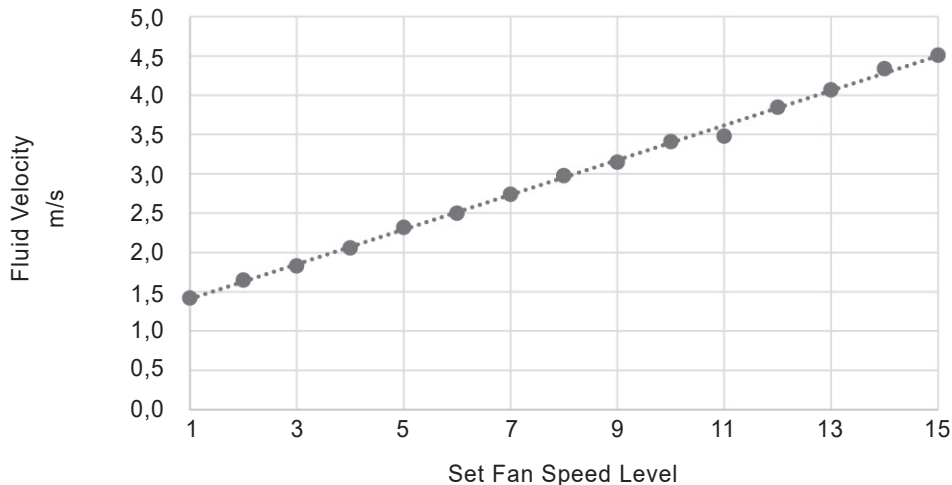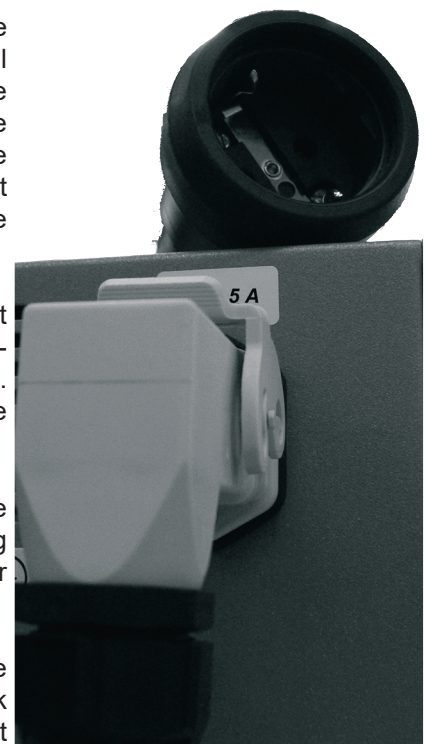
### 3.2.6 Connecting a Cool Line Refrigerator (2.00.321 C model version only)

The custom-designed power socket at the back of the enclosure (see the photograph to the right), which is designed for connecting a Cool Line series refrigerator to the unit, can be enabled or disabled with the “Cool Line” option of Configuration Menu 2. When it is disabled, the refrigerator is switched off and there is a “Cool Line disabled” message on the standard display, alternating with the status message “Exhaust Air OK” on the first line. Acceptable specifications of a refrigerator are 230 V AC, 50 Hz, max. 5 A.

 **Important** The 2.00.321 C model version of the exhaust air control unit ships with a dedicated adapter cable matched to the custom-designed power socket (see the photographs to the right). Power may be supplied to the cabinet **exclusively** with the factory-installed socket and shipped, dedicated cable.

 **Important** As soon as an error condition is detected, power to the power socket is switched off, which is equivalent to turning the refrigerator off. The same holds when the exhaust air monitoring is disabled in the Configuration Menu 2.

 **Info** When enabled, the special power socket at the back of the exhaust air control unit checks whether it receives a loop-back signal from the connected storage cabinet. If it is enabled, but does not detect a loop-back signal, a message “Error Cool Line” is displayed at the second line of the screen.

### 3.2.7 Enabling an Audible Alarm in Case of Extraction Airflow Error

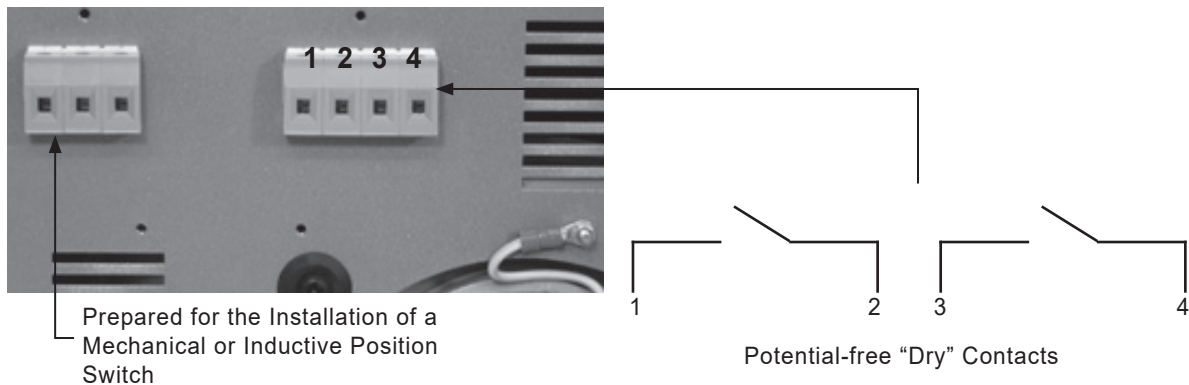
When the “Airflow Signal” submenu was selected, the ↑, ↓ cursor keys are used to select the desired behavior of the audible alarm signal that sounds with an extraction airflow error indicated on the screen. The selected item is saved and applied when you press START/SET.

The “No Delay” item means that the alarm will sound immediately when an error message appears on the screen. By selecting “Off,” you disable the alarm sound altogether. The time items (30 sec, 1 min,


1.5 min, 2 min, 3 min), when selected, will trigger the sound with a delay of 30 sec or 1 min etc. from the moment the error message is displayed.

### 3.3 Potential-free (“Dry”) Contacts

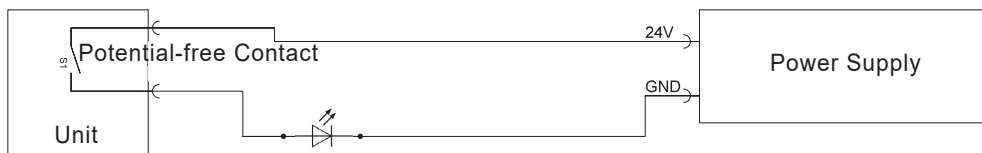An error condition causes two normally open, potential-free contacts to be closed; for individually locating errors, their closure can be recorded at the switching panel from the backside of the enclosure (see above).

Confirming the error and its correction by pressing the START/SET key (see Sect. 3.1) causes the two contacts to be reset (i.e., opened).

 **Important** The potential-free circuit (only one contact reproduced in the diagram) is designed for **max. 24 V, 2 A**. If these values were exceeded, the unit may be irrevocably destroyed. Additionally, any such damage would not be covered by the product warranty.

### 3.4 Setting the Pressure Difference

At a metering orifice, the suction line from the storage cabinet is equipped with a pitot probe. The pressure difference it senses is measured by a differential pressure meter via the two transparent tubes routed to it. It rises as the airflow through the suction-side extraction line on site increases. With restricted or blocked airflow, the pressure difference sensed by the probe drops. If it drops below the threshold set at the differential pressure meter, a signal that an error condition has developed is sent into the enclosure for further processing via the cable in the blue protective conduit. For technical reasons (hysteresis), the threshold from which the monitor cancels the error condition when the pressure difference rises again, is a little higher than the threshold that triggers an error. Therefore, make sure at all times that the settings for the pressure difference and the fan speed of the air extraction system on site are matched to each other in such a way that both an error is indicated when a restriction is detected, and the error is canceled by the unit when the restriction is cleared. There is no upper threshold for a maximum pressure difference.

The value of the pressure difference threshold can be set at the meter from 20 Pa to 300 Pa. This is done by simply removing the transparent cover and using a screwdriver in the provided slot to turn the selector. The factory-set default threshold of the pressure difference is 70 Pa.

Setting the pressure difference threshold for monitoring the airflow requires the following steps:

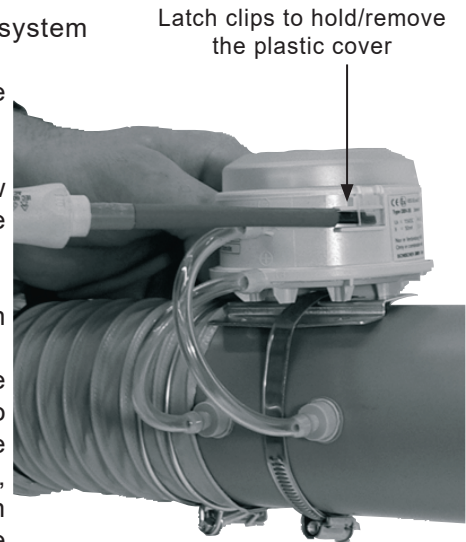
1. Connect the exhaust air monitor properly to the extraction system (outlet nipple) on site.
2. Set the fan speed to the desired extraction airflow (see Sect. 3.2.5).


Caution

In a cabinet for flammable liquids, the extraction airflow has to be sufficient to replace the interior volume of the cabinet 10 times every hour.

3. Loosen the plastic cover of the differential pressure meter on the suction hose (refer to the photograph to the right).
4. Using an appropriate screwdriver in the turning selector at the center of the differential pressure meter, set the threshold to make "Exhaust OK" appear on the screen (reduce pressure difference). If no error is displayed on the screen at the start, the pressure difference setting may first be increased until an error is displayed, then slowly reduced again. The higher the pressure difference threshold is set, the quicker / more sensitive the meter will respond by throwing an error message.
5. Using appropriate means on site (e.g. by covering the exhaust outlet with your hand), interrupt the extraction airflow temporarily.
6. If set correctly, the unit will trigger the audible alarm and display the error message on the screen. After the airflow interruption is cleared, the audible alarm should stop and the "Confirm error" message should appear on the screen.
7. Reinstall the plastic cover on the differential pressure meter.
8. Confirm the error message on the screen by pressing the START/SET key to re-open the dry contacts.

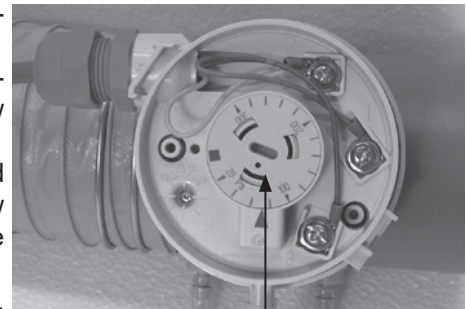Latch clips to hold/remove the plastic cover

Selector for setting the pressure difference threshold

### 3.5 Operation on Battery, Switching on and off with and without Battery

In the event of power supply failure, the control unit can be operated on the internal battery (accumulator) for approx. 3 hours.

This is why, provided the battery is sufficiently charged, pressing the On/Off switch will not turn the unit off. When the switch was pressed, the battery will continue to supply power to it. To turn the monitor off definitely, i.e. to disconnect power to it from both the mains network and the battery, hold down the STOP/RESET key for approx. 5 seconds when the standard display shows on the screen. This causes the message "Please disconnect power to unit!" to appear. Do as you are asked to: switch the monitor off using the On/Off switch and unplug the mains plug. Wait a short time for the monitor to switch itself off.

If the unit is switched off without holding down the STOP/RESET key first, this counts as a power supply failure. It will trigger an error message and close the dry contacts (see Sect. 3.1 and 3.3). In case the On/Off switch was pressed first, the monitor will indeed switch itself off when the STOP/RESET key is then held down. However, the error message that this will cause has to be confirmed when it is switched on again.

If the battery is charged to a sufficient level. it is possible to switch the unit on while it is not connected to mains power, by pressing the START/SET key. If the battery voltage drops below a factory-set threshold, the monitor switches itself off automatically. Turning on and operating the module without mains power will trigger a power supply error message as well, which will have to be confirmed.

### 3.6 Online Monitoring over the Internet with DÜPERTHAL connect

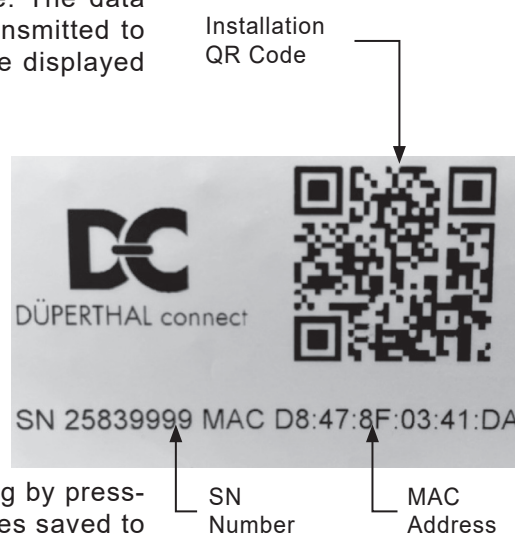
Using DÜPERTHAL connect, you can follow online the operating status and any malfunctions or alerts of the monitoring and air extraction units. A connection to DÜPERTHAL connect is possible only with units that are capable of radio communication (P/N 2.00.321 C, 2.00.322).

Furthermore, a Bridge is required as the communication gateway in order to connect the unit to DÜPERTHAL connect. When a Bridge is within reach, the unit can be registered with DÜPERTHAL connect immediately. To do so, perform the following steps:

1. On the DÜPERTHAL connect home page, access the “Monitoring” menu item.
2. Use the “Add monitoring” button to start the installation process.
3. As the first step, select the cabinet in question.
4. Then, select the appropriate hub in the selection listing or add a new one with the help of the “Create new hub” button.
5. A new hub can be added by scanning the QR code present on the name plate tag of the unit itself or by entering its SN (serial number) and MAC address indicated on the name plate manually.
6. Select the desired functionality of the installation site. The data pertaining to the selected functionality will then be transmitted to the DÜPERTHAL connect online monitoring and will be displayed at the installation site.

If you enter the SN and the MAC address of the unit manually in DÜPERTHAL connect, separate both by a semicolon “;” (in the case of the example in the margin, the required input is: 25839999;D8:47:8F:03:41:DA).

### 3.7 Error and Menu Change Reporting

Error conditions and changes applied to the configuration menus are recorded and saved to memory. They can be viewed after pressing the ⇐ arrow key in the standard display (see Display 11). You can scroll through the displayed listing by pressing the ↑ and ↓ arrow keys. The maximum number of entries saved to memory and displayed in this listing in chronological order is 50. When the maximum number is reached, any further report will overwrite the oldest one in memory. Deleting the entries in memory is possible from the Supervisor Menu.

A report entry contains information on the type of entry, i.e. error or menu change, the date, and the time of day. The title line of the listing shows the number of the report currently looked at (before the slash) and the total number of entries in memory (after the slash). If the entry is an error report, the error code (E1 or E2) is shown. These error codes have the following meaning:

- E1 Power Supply Error
- E2 Exhaust Air Error

The letter “M” stands for the report of a modification of the configuration menus. As is indicated by the symbolized arrow to the right, pressing the ⇐ arrow key, when this entry is selected, will show additional information. This information is a list of all settings of the configuration menus. All the options are given with the indication of their status, changed or unchanged at the moment the modification record was generated. The record for a modification is generated only when the operator quits the configuration menus, and it includes all settings, whether changed at that point in time or remained as is.

* Report 018/42 *
E1 22.06.08 13:32
>M 22.06.08 13:41->
E2 24.06.08 08:09

Display 11

The following overview is meant to assist you in interpreting the change reports.

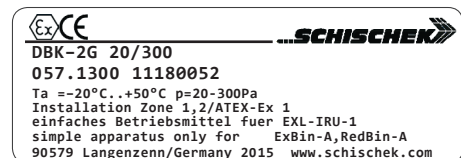
- Date/Time <> Date/time before do not equal after change
- Language = Set language before equals after change
- Exh Monitor 1->0 The exhaust monitoring setting was set from “on” (1) to “off” (2)
- Speed 12->9 Fan speed was changed from level 12 to level 9
- Cool Line 1->1 The setting for the refrigerator socket was left at “on” (with 2.00.321 C model only)

The report screen is quit by pressing the STOP/RESET key, or it returns automatically to the standard display after a short time of inactivity.

#### 4 Technical Specifications

- Power Supply Voltage 230 V
- Frequency 50 Hz
- Current Rating 0.6 A
- Motor Power 100 W
- Revolutions 2810 /min
- Fluid Velocity max. 4.51 m/s
- Volume Flow Rate max. 63.5 m<sup>3</sup>/h
- Weight approx. 9.0 kg
- Dimensions (enclosure edges) 260 x 310 x 275 mm
- Ingress Protection IP20
- ATEX Rating EX II (3) G [c II B T4]

see also the component name plate, as to the right

#### 5 Troubleshooting

In the course of using the appliance, malfunctions may occur. The table below is meant to help troubleshoot and correct malfunctions. If a malfunction cannot be traced to one of the possible causes or the correction given does not solve the issue, then use of the appliance has to be discontinued. Consult with the distributor on troubleshooting an repair.

Malfunction	Possible Cause	Verification of Cause	Correction of Malfunction
Appliance does not run	On/Off Switch Power Supply Cord Overvoltage Switch Power Socket at the Back (Cool-Line cabinet)	<ul style="list-style-type: none"> <li>· Is the appliance switched on at the On/Off switch?</li> <li>· Is the power supply cord damaged?</li> <li>· Is the power supply cord properly connected?</li> <li>· Are the fuses inside the appliance o.k.?</li> <li>· Is a device connected that may consume current at more than 5 A?</li> </ul>	<ul style="list-style-type: none"> <li>· Switch the appliance on at the On/Off switch.</li> <li>· Replace the supply cord with a new one.</li> <li>· Connect the supply cord to the intended socket.</li> <li>· Replace the fuses (2 x 1.25 A slow blow).</li> <li>· Disconnect this device from the socket and check the fuses and replace them as needed (2 x 5 A slow blow).</li> </ul>
Fan runs, but display screen does not work	Display Electrics/ Electronics	· Malfunction cannot be corrected by user.	· Consult with the manufacturer/distributor on inspection/repair.
Display screen works, but fan does not run	Fan Speed Setting	· Is the fan speed set to level 0?	· Set the fan speed to a higher level (see Sect. 3.2.5).
Power Supply Error (E1) shows on screen	<i>Possible Cause, Verification, and Correction as under “Appliance does not run” above.</i>		

Malfunction	Possible Cause	Verification of Cause	Correction of Malfunction
Exhaust Airflow Error (E2) shows on screen	Inlet or Outlet Hose  Differential Pressure Meter/Switch  Switching Amplifier	<ul style="list-style-type: none"> <li>Is the suction or the extraction hose of the appliance blocked (airflow restriction) or very dirty?</li> <li>Are the measuring tubes connecting the sleeve to the meter/switch properly connected?</li> <li>Are the measuring tubes connecting the sleeve to the meter/switch damaged?</li> <li>Are the measuring tubes connecting the sleeve to the meter/switch connected to the correct nipples?</li> <li>Is the differential pressure meter/switch properly set?</li> <li>Is the spiral hose connected tightly to the sleeve?</li> <li>Is the differential pressure meter/switch wired properly?</li> <li>Is a clapping noise heard when the differential pressure switch cuts in or out?</li> </ul>	<ul style="list-style-type: none"> <li>Clean the hose and/or clear the restriction.</li> <li>Connect the measuring and connecting tubes to both the meter/switch and the sleeve in the extraction line.</li> <li>Replace the measuring and connecting tubes with new ones.</li> <li>Connect the tubes as intended: nipple closer to the cabinet of the sleeve to P1 (+) of the meter/switch, nipple closer to the appliance of the sleeve to P2 (-) of the meter/switch.</li> <li>Open the differential pressure meter/switch and set it properly (see Sect. 3.4).</li> <li>Tightly set the spiral hose onto the sleeve using the clamp.</li> <li>Open the differential pressure meter/switch and wire it correctly: blue wire to 1, yellow to 2, red to 3.</li> <li>Consult with the manufacturer/distributor on inspection/repair.</li> </ul>
Despite irregular extraction airflow no airflow error (E2) on screen	Configuration Menu 2 Differ. Pressure Meter/Switch	<ul style="list-style-type: none"> <li>Is the feature for monitoring the extraction airflow turned off?</li> <li>Is the differential pressure meter/switch properly set?</li> </ul>	<ul style="list-style-type: none"> <li>Set the menu option for the exhaust airflow monitor to "on".</li> <li>Open the differential pressure meter/switch and set it properly (see Sect. 3.4).</li> </ul>

## 6 Contact Details

DÜPERTHAL Sicherheitstechnik GmbH & Co. KG  
 Frankenstr. 3  
 63791 Karlstein, Germany  
 Web: [www.dueperthal.com](http://www.dueperthal.com)

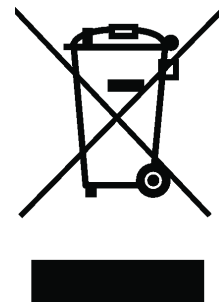
Tel.: +49 6188 9139 0  
 Fax: +49 6188 9139 121  
 E-mail: [info@dueperthal.com](mailto:info@dueperthal.com)

We reserve the right to change technical specs of the unit without prior notice.

Pursuant to the directive 2012/19/EU on Waste Electrical and Electronic Equipment (so-called WEEE Directive), equipment that was manufactured or distributed by ourselves may be returned to us. To discuss the exact procedure we will follow, please contact us with the details above.

We also declare that the equipment manufacture conforms to the directive 2011/65/EU on the restriction of the use of certain hazardous substances in electrical and electronic equipment (so-called RoHS Directive).

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